



Genemed Synthesis, Inc.  
6203 Wood Lake Center Dr., Bldg. 2, San Antonio, TX 78244, USA  
**Toll free** (800) 344-5337; **Phone:** (210) 745-5988; **Fax** (210) 745-5992  
**Email:** [info@genemedsyn.com](mailto:info@genemedsyn.com), **Website:** [www.genemedsyn.com](http://www.genemedsyn.com)

Product data sheet

---

### Blocking Buffer Concentrate

**Cat. No.** EW-80060

**Size:** 1 x 50 ml

**Form:** 10X Solution

**Storage:** Store at 4<sup>0</sup>C

---

#### Introduction

Antigen or antibody coating on ELISA plates is commonly used as the first step in ELISA. This is followed by the blocking of the excess protein binding sites in the microwell not occupied by the coated antigen/ antibody. Genemed synthesis Inc's Blocking Buffer has been specially formulated to serve two important functions: to block the excess protein binding sites on ELISA plates and to stabilize the coated antigen or antibody on the plate. The ELISA plates blocked with this solution can be stored dried at 4<sup>0</sup>C for an extended period of time.

#### Form

The blocking buffer concentrate is formulated using BSA and contains 0.02% merthiolate as preservative. **It should not be used in applications where BSA interferes with the assay.** (gelatin containing blocking buffer, catalog no. EW-80061, or milk-based buffers are also available from Genemed Synthesis)

#### Recommended Use

Dilute 1:10 with distilled water before use. Add 150-200  $\mu$ l of diluted solution per well and incubate plates for a minimum of 2 h at 37<sup>0</sup>C or 3-24 h at room temperature. After incubation, aspirate solution and air dry ELISA plates without washing. After drying, plates can be stored at 4<sup>0</sup>C. For some applications, it may be necessary to block at 1:5 dilution to reduce background.

The blocking buffer is most compatible with high protein binding ELISA plates (Cat. No. EW-80011). It is not recommended for other applications such as membrane blocking (Western blotting, etc.).

**MSDS:** A Material Safety Data Sheet is not required for this product. The product does not contain any hazardous components above 1% or any carcinogens above 0.1% as defined in 29 CFR 1910.1200, the OSHA Hazard Communication Standard. The product contains natural purified proteins in a buffer containing 0.1% or less sodium azide. MSDS can be consulted for these items. Please note that although the product is defined as not hazardous, it is still advisable to follow prudent laboratory practices when handling laboratory reagents.

This material is sold for research purposes only and is not required to appear on the TSCA inventory. It is not intended for food, diagnostic, drug, household, agricultural or cosmetic use. Its use must be supervised by a technically qualified individual experienced in handling potentially hazardous chemicals.

*This product is for in vitro research use only.*

---

#### Related items available from GSI

Catalog#	ProdDescription
EW-80012	High binding ELISA Strips plates (8 wellsx12 strips)
EW-80050	ELISA Plate Coating buffer concentrate (10X)
EW-80060	ELISA Plate Blocking Buffer concentrate (10X) BEW-based
EW-80062	ELISA Plate Blocking Buffer concentrate (10X) milk-based
EW-80070	Antibody and Conjugate Diluent for ELISA concentrate (10X)
EW-80080	Wash buffer concentrate (20X) for ELISA
EW-80081	Phosphate Buffered Saline (PBS, pH 7.4) concentrate (20X)
EW-80091	TMB substrate (1-component) for ELISA
EW-80100	Stop solution for TMB substrate (ELISA) concentrate (10X)
EW-80110	Single solution TMB substrate (membrane) solution for western blot
EK-80160	Rabbit Serum Antibody detection ELISA kit, Qualitative (mouse, goat, sheep, monkey, human, chicken kits are also available)

EW-80060

Rev. 80609F